

Transfer Data from CI-340 to PC:

1. Connect the CI-340 to the computer and power it on.
2. Create a driver for the CI-340, using the inf-wizard.exe. This is located in the **libusb-win32 bin 1.2.6.0** folder on the software cd or can be downloaded at <http://sourceforge.net/projects/libusb-win32/files/libusb-win32-releases/1.2.6.0/libusb-win32-bin-1.2.6.0.zip/download>
 - a. Double-click on the inf-wizard.exe to start the driver creation wizard.
 - b. Click “run” to run the application even though the publisher could not be verified.
 - c. Now, in the inf-wizard application, click “Next” after making sure the CI-340 is connected to the computer and powered on.
 - d. Select the correct device from the list. The Vendor ID is 0x0425 and the product ID is 0x0500. The device description may say “CI-340”. Select the 340 from the list and click “Next.”
 - e. Change the Device name to CI-340 if it is not already so. Click “Next” after configuring the device.
 - f. You will be prompted to save the CI-340.inf file. Remember where you save the file-this is the driver for the CI-340.
 - g. After saving the file, in the inf-wizard, click the “Install Now” button to install the driver you just created.
 - h. Click “Yes” to allow an unknown publisher to make changes on the computer.
 - i. Click “Install this driver software anyway” when windows can’t verify the publisher.
 - j. The installation process can take a few minutes to complete. When it is successful, click “Okay.”
3. Now, install and open the data transfer software by double-clicking the C340dfv6010.exe. The application to download data from the CI-340 can be found on the web at <http://cid-inc.com/support/ci-340-support/340software>
 - a. If the publisher cannot be verified, click to Run the application anyway.
 - b. In the software application “CID Download Data from CI-340”, click “File” and “Open” to see the files on the CI-340.
 - c. Select the files from the list to transfer to the computer and click “Open.”
 - d. Save the file to your computer and re-open it with a spreadsheet program such as Microsoft Excel. The file type will be .csv or comma separated value.

With **Windows 8**, the driver signature enforcement may need to be disabled to connect the CI-340 to the PC. For instructions, please visit: <http://www.howtogeek.com/167723/how-to-disable-driver-signature-verification-on-64-bit-windows-8.1-so-that-you-can-install-unsigned-drivers/>

