

Replacing Previous Versions of the CI-600 Driver:

If this system already has a previous version of CI-600 driver installed these steps must be followed when installing the new version of software. After downloading or installing the software (www.cid-inc.com/Software/CI-600):

1. Plug-in the CI-600 instrument.
2. Click the **Start** button.
3. Right click **Computer** and select **Properties**.
4. Click on the **Device Manager** link.
5. Expand the **Imaging Devices** and right click on **CanoScan LiDE 30/N1240U**.
6. Click on **Update Driver Software...**
7. Select “**Browse my computer for driver software**”.
8. Select “**Let me pick from a list of device drivers on my computer**”.
9. Click the **Have Disk** button.
10. Type (or browse to select) the path of the CI-600 driver software.

By default the path is...

C:\Users\YourUserName\Documents\CI-600 Driver

11. Select **CI-600 Root Scanner** from the list of drivers.
12. Click **Next**.
13. If dialogs appear warning that this driver has not been signed by Microsoft, click **Continue** or **OK** to install the driver.
14. If dialogs appear asking if you want to replace any files, click **OK** , **Replace** or **Yes** to continue.
15. Reboot the computer.
16. Consult the CI-600 Instruction Manual for calibration and scanning instructions.